

# Determination of Silicon Carbide in Recycled Catalysts



Silicon Carbide filter

This semi-quantitative method for determining silicon carbide (SiC) content in mixed batches of recycled emission control catalysts uses thermogravimetric analysis and total carbon analysis.

#### **Determination of %SiC**

SiC is most often found in Diesel Particulate Filters (DPF's) that are part of an emissions catalyst system. This method to determine the %SiC is based on a published standard (Ref. No. ISO 21068-2:2008) and employs:

- LECO 701 Thermogravimetric Analyzer (TGA) for automated Loss on Drying/Ignition determination
- LECO CS600 Carbon/Sulfur Analyzer for coulometric determination of total carbon

#### Perform Analytical Analysis

- Run LOD (Loss on Drying) @250°C
- Run LOI (Loss on Ignition) @650°C
- %C Volatile (free carbon) is LOI LOD
- Use LECO C analyzer for %C Total


## $\rightarrow$

#### Calculate %SiC

- %C in SiC = %C Total %C Volatile
- %SiC = (%C in SiC) ÷ 0.2995

#### **Method Validation**

There is excellent correlation between this method and research grade X-ray diffraction methods.

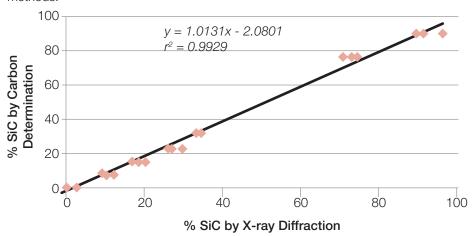

#### **About Us**

BASF's Catalysts division is the world's leading supplier of environmental and process catalysts. The group offers exceptional expertise in the development of technologies that protect the air we breathe, produce the fuels that power our world and ensure efficient production of a wide variety of chemicals, plastics and other products, including advanced battery materials. By leveraging our industry-leading R&D platforms, passion for innovation and deep knowledge of precious and base metals, BASF's Catalysts division develops unique, proprietary solutions that drive customer success.

#### BASF - We create chemistry

#### Contact us today

## BASF Catalysts Headquarters

BASF Corporation, Catalysts Division 25 Middlesex / Essex Turnpike Iselin, New Jersey, 08830, USA

Tel: +1-732-205-6060

#### Americas

Tracy Ware-Justice (973) 715-5677 tracy.ware-justice@basf.com

Keith Allison (732) 762-4043 keith.allison@basf.com

#### Asia-Pacific

David Zhao +86 188 0190 2818 david.zhao@basf.com

### Europe, Middle East, Africa

Bob Smallwood +44 776 924 8526 bob.smallwood@basf.com

Dave Foley +44 7831 806094 david.foley@basf.com

Stephen Dew Nash +44 7775 221957 stephen.dew-nash@basf.com

Although all statements and information in this publication are believed to be accurate and reliable, they are presented gratis and for guidance only, and risks and liability for results obtained by use of the products or application of the suggestions described are assumed by the user. NO WARRANTIES OF ANY KIND, EITHER EXPRESS OR IMPLIED, INCLUDING WARRANTIES OF MERCHANTABILITY OR FITNESS FOR A PARTICULAR PURPOSE, ARE MADE REGARDING PRODUCTS DESCRIBED OR DESIGNS, DATA OR INFORMATION SET FORTH. Statements or suggestions concerning possible use of the products are made without representation or warranty that any such use is free of patent infringement and are not recommendations to infringe any patent. The user should not assume that toxicity data and safety measures are indicated or that other measures may not be required.

© 2017 BASF

BF-10564 09/17